



Fast G.C. Side Cover & Cryo Mixer Sub.

Subassembly

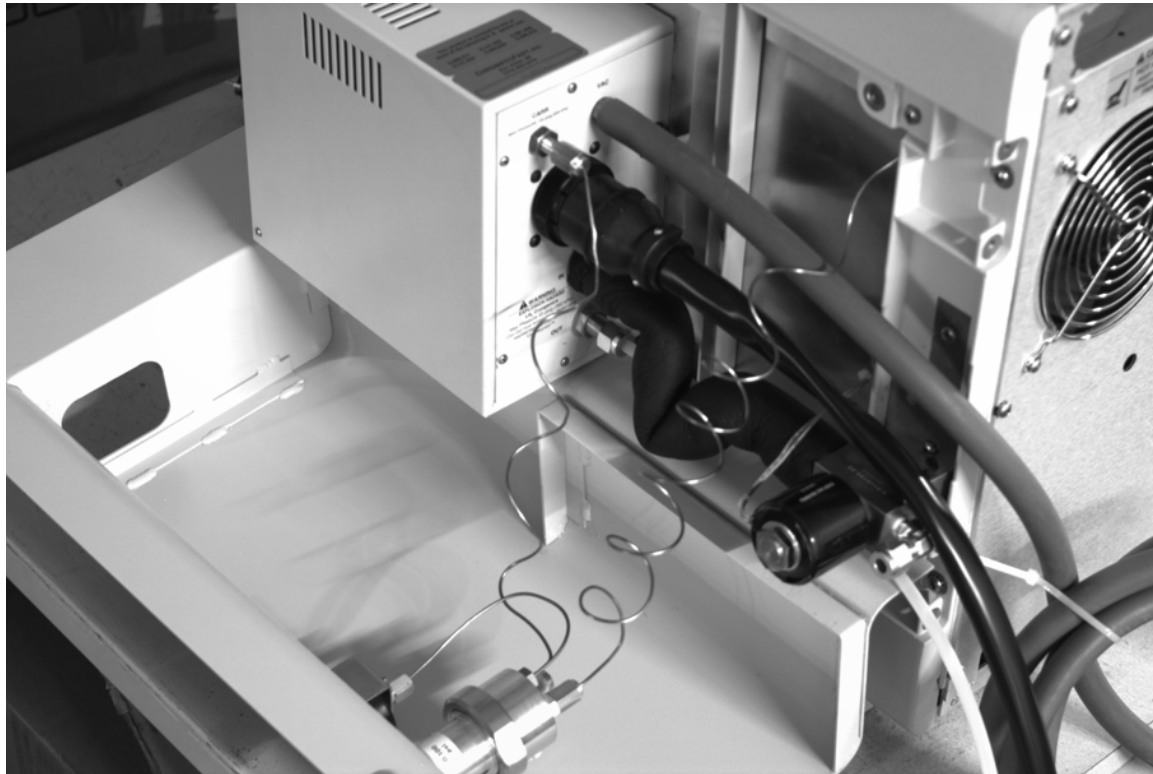

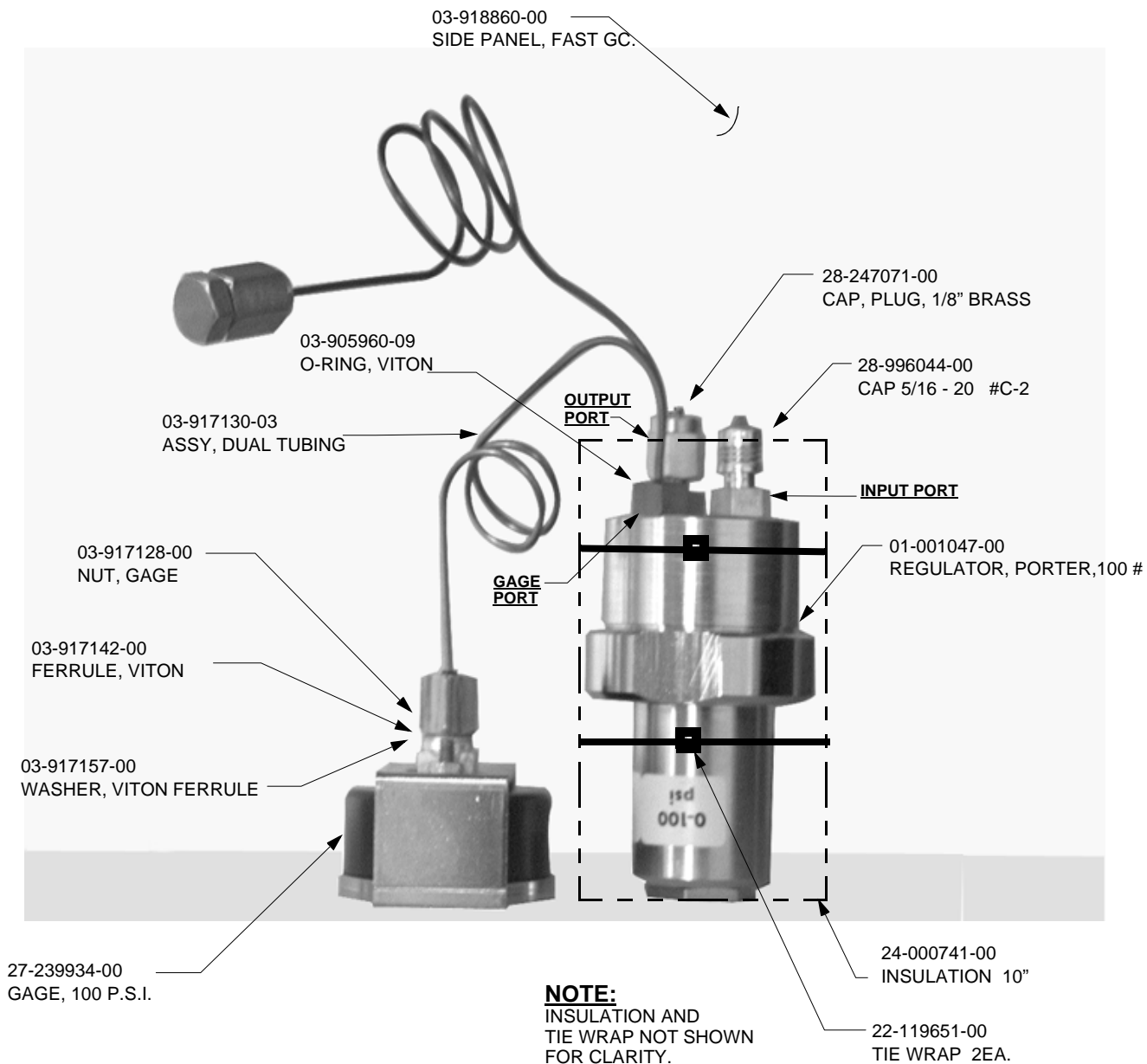

NOTE:
USE T-4343 FOR LABEL PLACEMENT

88-494510-00
PIPE INSULATION (LOCATION ONLY,
INSULATION NOT SHOWN. INSUL. TO
COVER ALL FITTINGS)

28-691946-00
1/4" BRASS TEE


28-995813-00
VINYL CAP .625 DIA.

28-995806-00
VINYL CAP, .125 DIA.

28-694540-01
REDUCING FERRULE,
1/4 NUT TO 1/8 TEE

03-925325-00
CRYO MIXER ASSY.

12-222060-10
SCREW, PHM 10-32 X 5/8 LG. SST

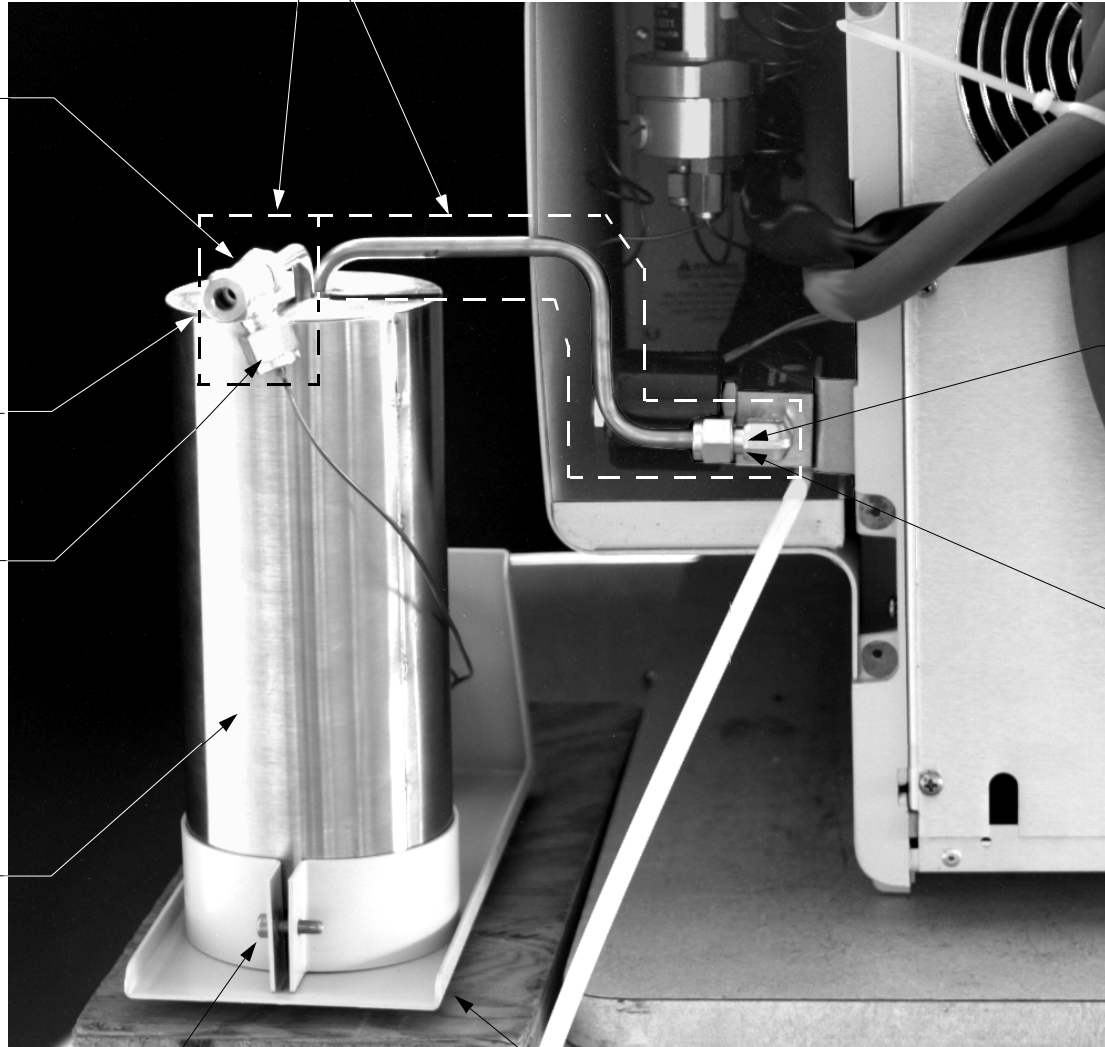

28-995808-00
VINYL CAP .250 DIA.

22-119650-00
TIE WRAP

23-608980-00
ADHEASIVE BACKED BUMPER 4EA

NOTES

1. FOR SHIPPING WITH FAST G.C. , DETACH MIXER FROM BASE & INSTALL CAP (28-995806-00) OVER TUBE END & PLACE BOTH PIECES IN SHIPPING BOX (INTEGRATOR), & RE-SEAL

03-925322-00
MIXER ASSY. BASE

SEE NOTE 2

88-189609-00
TFE TAPE, 1/4" WIDE

2X

28-849718-00
FITTING, 1/16 BARB TO
1/8 NPT, 90DEG.

27-239934-00
GAUGE, 100PSI

03-917128-00
NUT, 1/8"NPS - O-RING


03-917142-00
VITON FERRULE

03-917157-00
WASHER, FERRULE

03-947137-00
TUBE, GAGE ASSY.

NOTES

1. LOCATE FITTING SUCH THAT THE BARB PLUG (28-849973-00) POINTS DOWN.
2. POSITION REGULATOR SUCH THAT INLET PORT IS LOCATED ON THE RIGHT SIDE AS YOU LOOK AT THE GUAGE FACE AND THE REGULATOR KNOB.

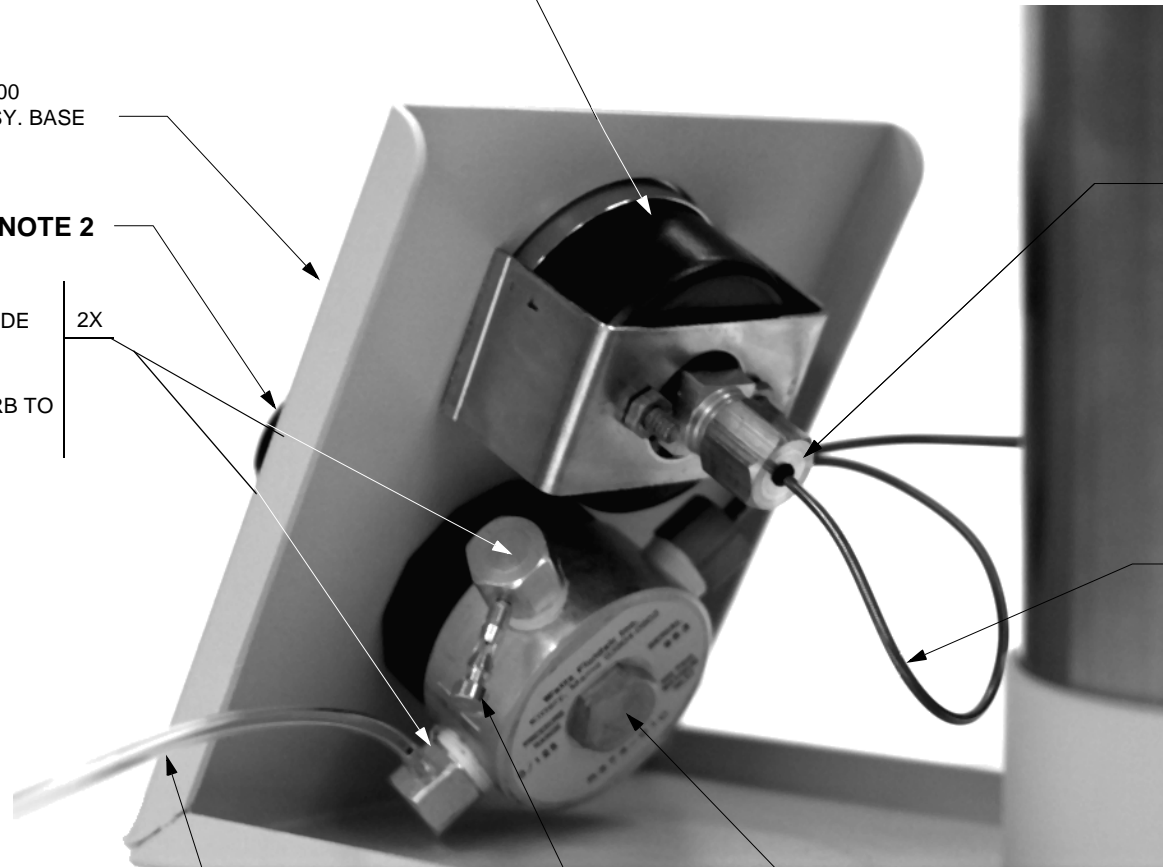

28-158923-00
TYGON TUBE 1/8' O.D.

27-189315-00
REGULATOR, 1/8", 2 WAY

28-849973-00
PLUG, BARB 1/16 W/POLY TUBE

SEE NOTE 1.