


Main Door Assembly

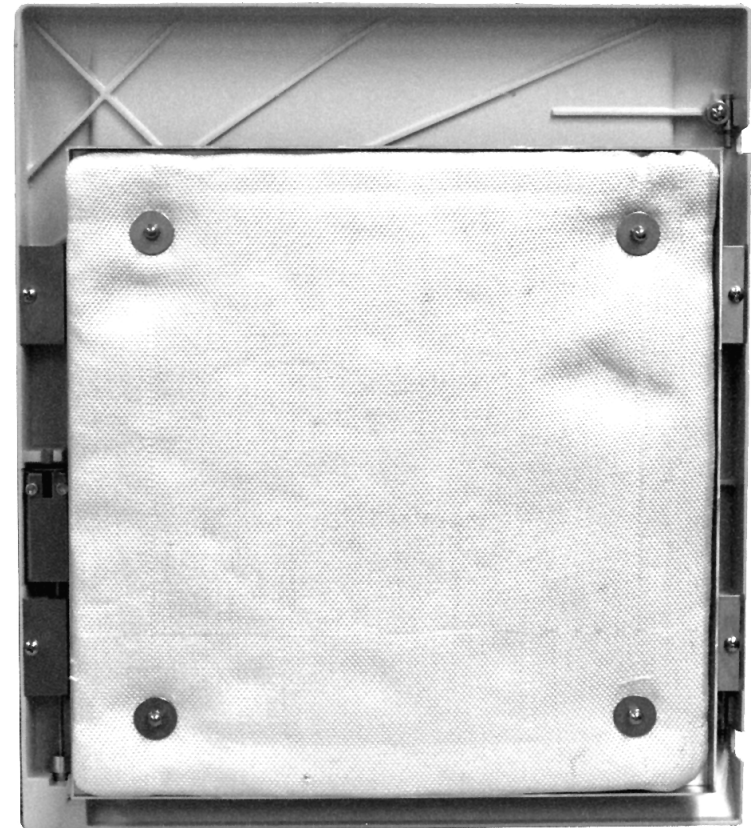

03-925208-01
MAIN DOOR

R

03-925311-01
PIN, HINGE, 2EA.

14-112007-00
WASHER, FLAT, 2EA.

12-901163-00
SCRW, 8-32 X 3/8, EXT.,LW, PH 2EA.

NOTES:

1. USE LOCTITE #222 ON BOTH SCREWS OF RETAINER LATCH. TIGHTEN SCREWS UNTIL SNUG OR LATCH PLATE MOVES. STOP IF PLATE MOVES.
2. ATTACH WIDE FLANGE OF OVEN DOOR SPRING ON LATCH SIDE.
3. FOR LATCH ROD ASSY. USE T# 4340

03-925283-01
PLATE, LATCH

03-925284-01
DOOR, RETAINER LATCH

03-925286-01
LATCH ROD

88-299284-00
LOCTITE #222
SEE NOTE 1

12-222008-08
SCRW, 8-32 X 1/2" SST, PH 2EA

17-699907-00
RIVET, 1/8 X 1/8, AL / AL 4EA.

NOTE:

TO ENSURE PROPER SEAL WHEN DOOR IS CLOSED, HANDLE DOOR ASSY. SUCH THAT NO DEPRESSION IS MADE IN INSULATION

03-925448-01
INNER COL. OVEN DOOR ASSY.

12-901132-00
SCREW, 8-32 X 3/8, EXT., PH, 4EA.

03-925287-01
SPRING, OVEN DOOR, 2EA.